

6.1 Parts Assembly

The following illustrations show most assembly parts and how they fit into the centrifuge.

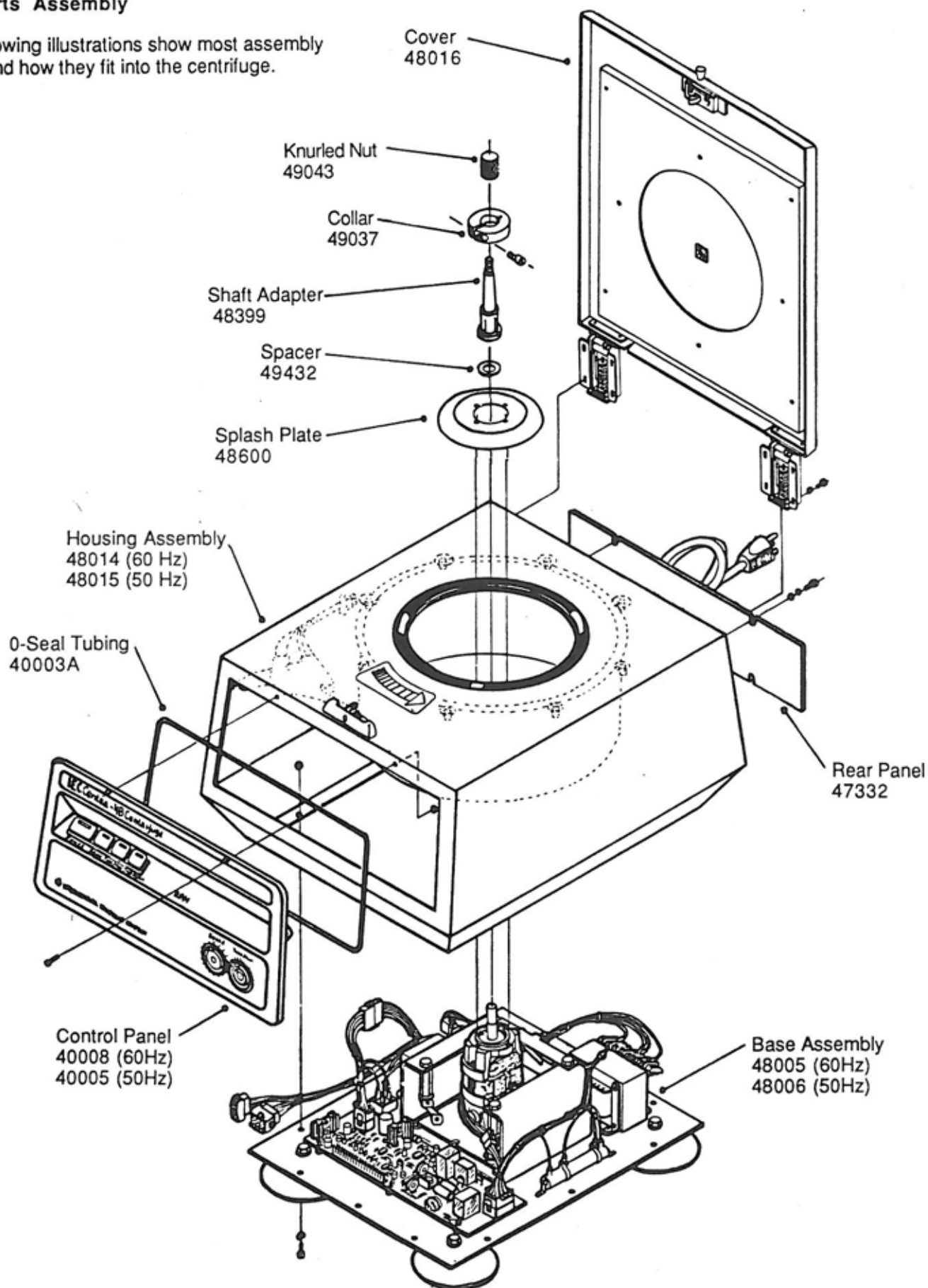


Figure 6-1 Centrifuge Assembly (1 of 3)

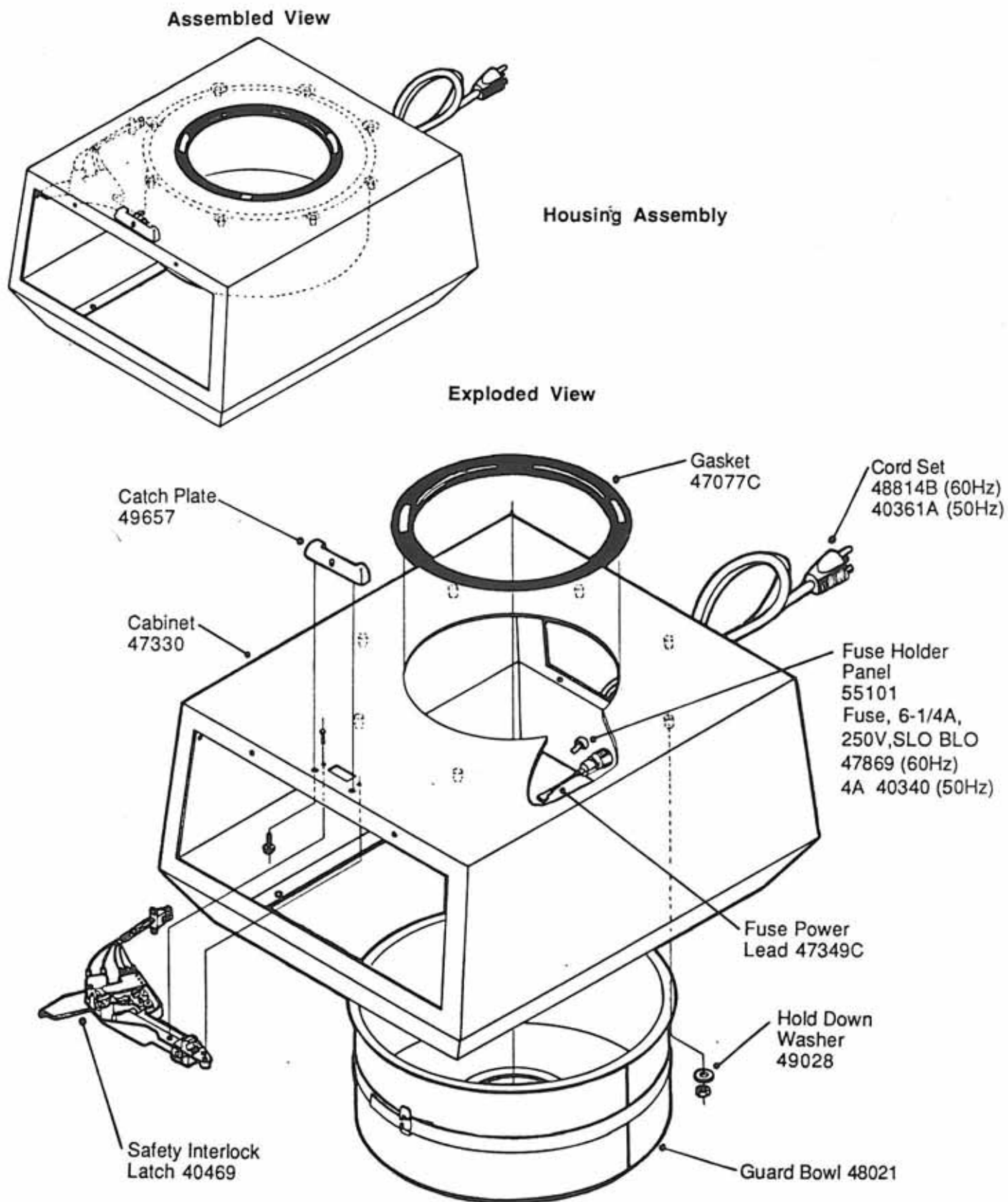


Figure 6-1A Centrifuge Assembly (2 of 3)

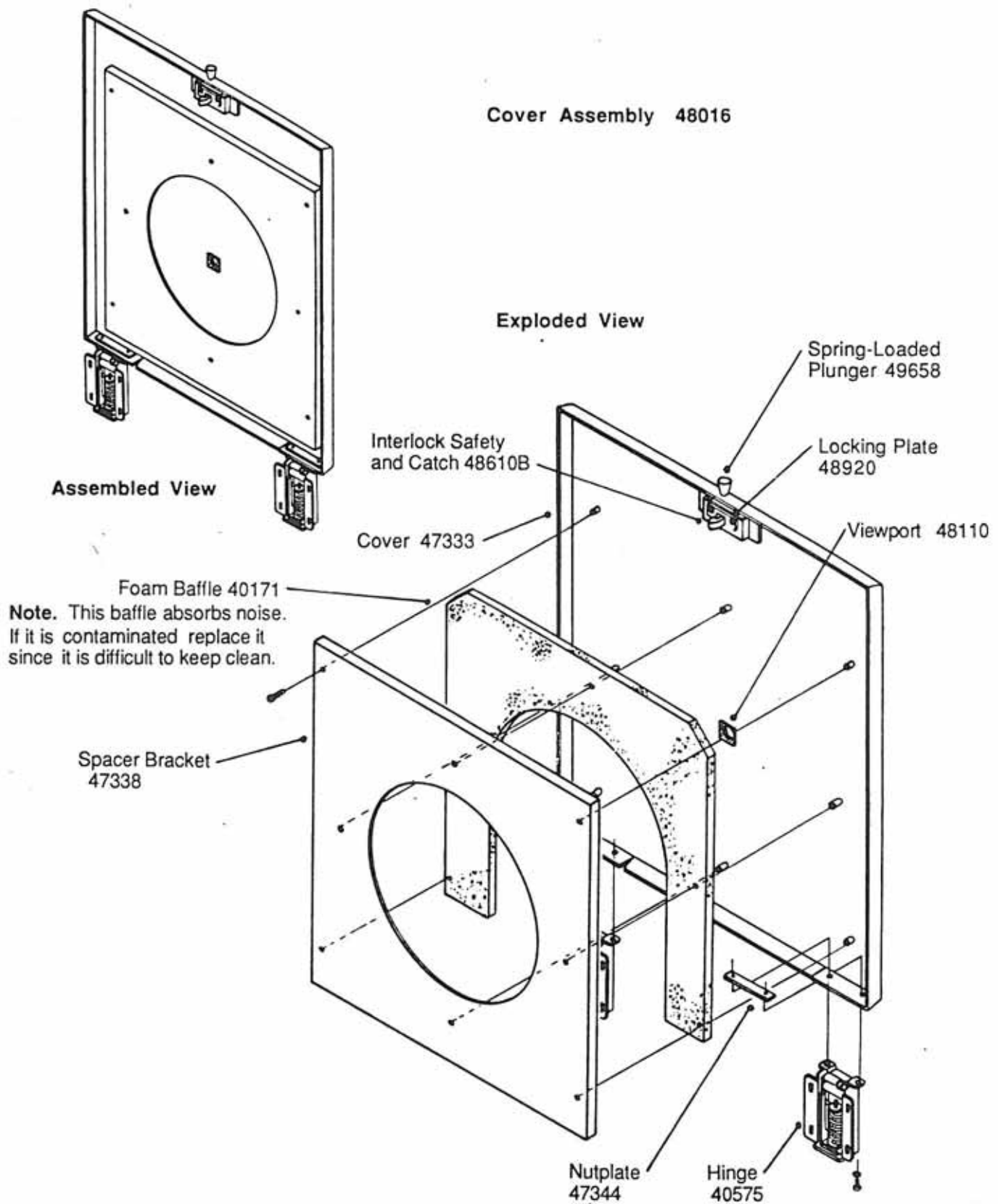


Figure 6-1B Centrifuge Assembly (3 of 3)

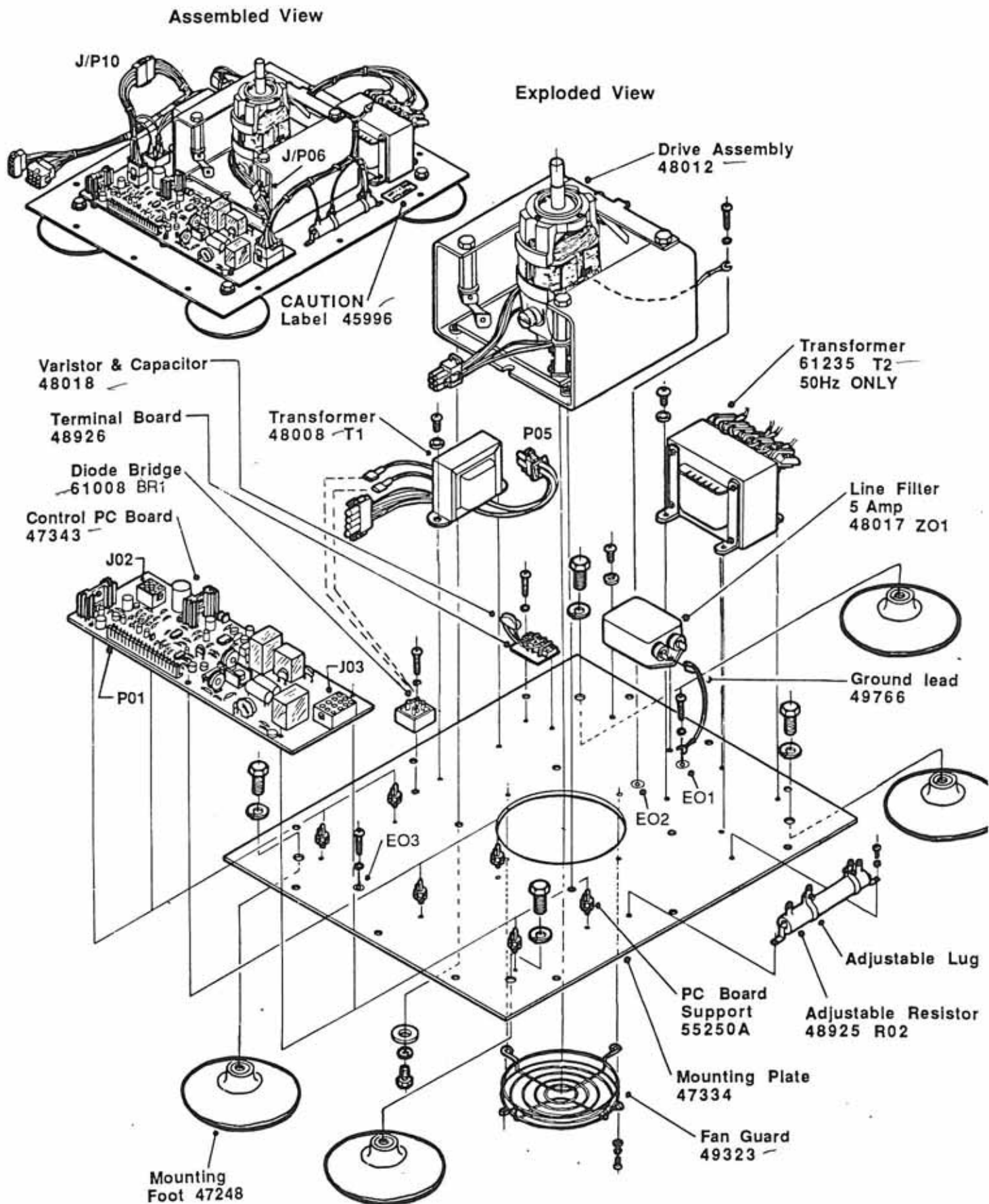


Figure 6-2 Base Assembly

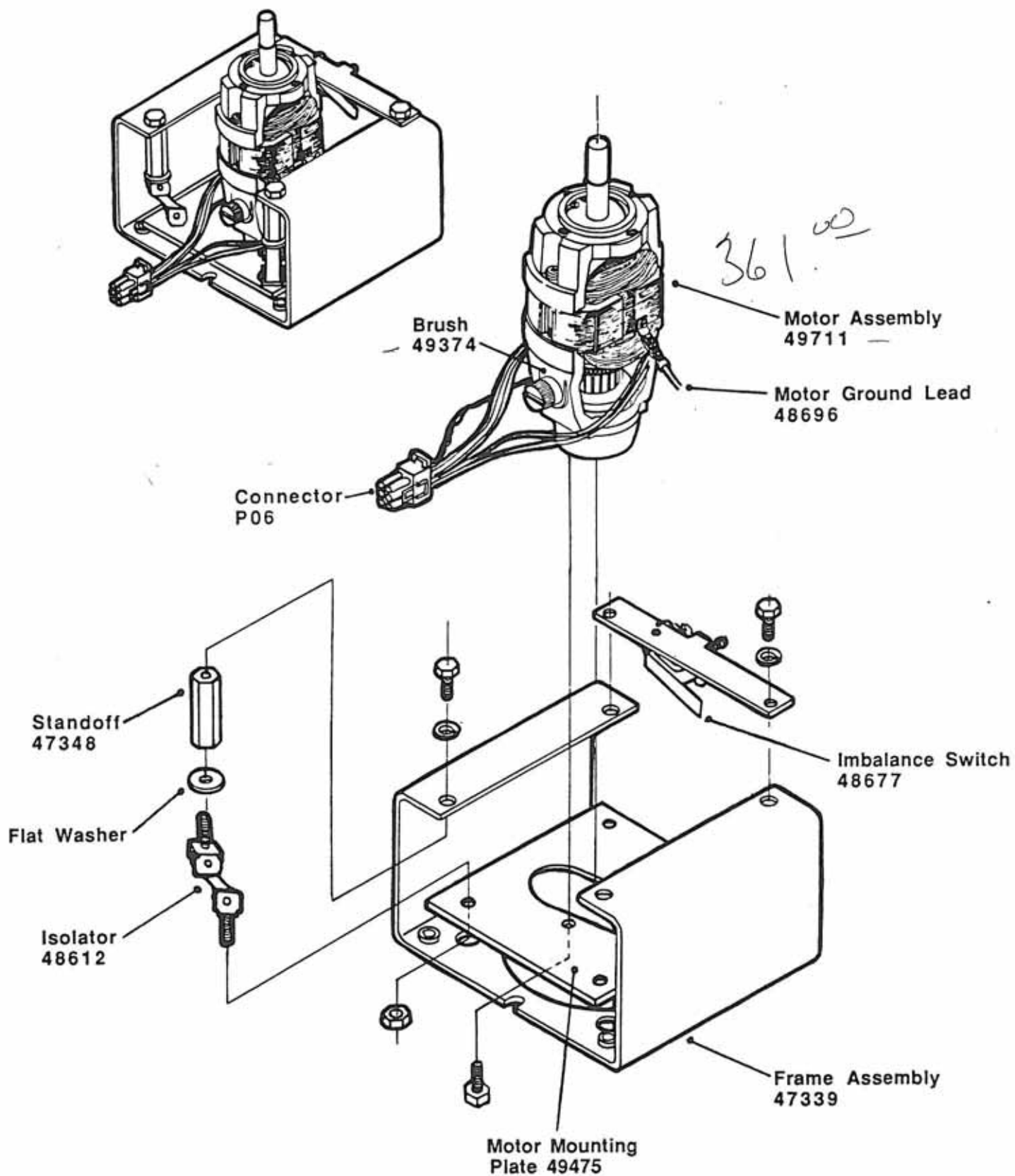


Figure 6-3 Drive Assembly 48012

6.2 Spare Parts

The following parts are recommended as spare parts.

Item	Description	Part No.	Quantity
1.	Brush Assembly	49374	One pair
2.	6-1/4 A, 250V fuse (for 100 or 120V,50 or 60 Hz units)	47869(Bussman MDA)	Two
3.	4 A , 250V fuse (for 200, 220 or 240V,50 or 60 Hz units)	40340(Littlefuse 326,004)	Two
4.	Foam Baffle	40171	One
5.	Rotor Locking Nut	49043	One