



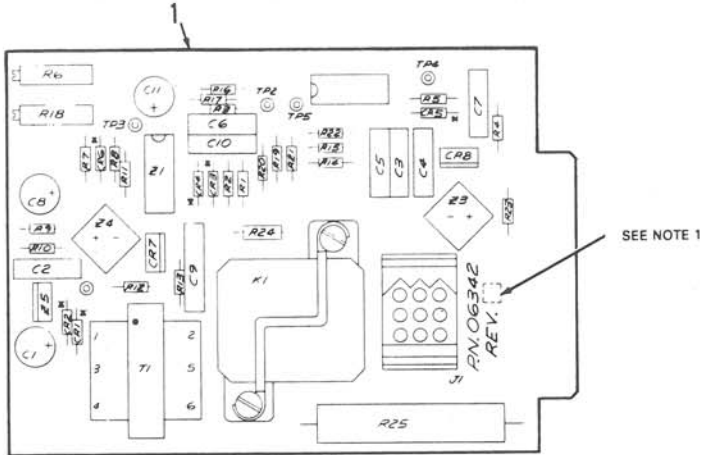

REF.	DWG.NO.	DESCRIPTION	REQ.
1	06320	PRINTED CIRCUIT BOARD ASSEMBLY	1
2	66141	CAPACITOR, SOLID ELECTROLYTE, DIPPED, 100 MF, 20V ± 10%	1
3	911006	CAPACITOR, CERAMIC, DISC, .1 MF, 50V	5
4	911002	CAPACITOR, CERAMIC, DISC, .01 MF, 1000V	1
5	911001	CAPACITOR, CERAMIC, DISC, .1 MF, 500V	1
6	923002	RECTIFIER, SILICON, 100V, IN4002	2
7	922501	DIODE, SIGNAL, SILICON, IN914	1
8	06334	CONNECTOR ASSEMBLY	1
9	65984	RELAY, 3 PDT, 23Vdc/115 AC, 10 AMP	1
10	900087	RESISTOR, COMP., CARBON 5%, 1/4W 4700 OHM	1
11	900069	RESISTOR, COMP., CARBON 5%, 1/4W 1000 OHM	4
12	900089	RESISTOR, COMP., CARBON 5%, 1/4W 5600 OHM	1
13	900063	RESISTOR, COMP., CARBON 5%, 1/4W 560 OHM	1
14	900063	RESISTOR, COMP., CARBON 5%, 1/4W 560 OHM	1
14	904007	RESISTOR, COMP., VARIABLE, CERMET, P.C. TYPE, 1000 OHM	1
15	900115	RESISTOR, COMP., CARBON 5%, 1/4W, 51K OHM	1
16	900115	RESISTOR, COMP., 5%, 1/4W, 2000 OHM	2
17	900083	RESISTOR, COMP., 5%, 1/4W, 3300 OHM	2
18	900047	RESISTOR, COMP., 5%, 1/4W, 150 OHM	2
19	900142	RESISTOR, COMP., 5%, 1/4W, 510K OHM	2

REF.	DWG.NO.	DESCRIPTION	REQ.
20	900016	RESISTOR, COMP., 5%, 1/4W, 10 OHM	1
21	900096	RESISTOR, COMP., 5%, 1/4W, 10K OHM	1
22	904013	RESISTOR, VARIABLE, CERMET, P.C. TYPE, 50K OHM	1
23	900086	RESISTOR, COMP., CARBON 5%, 1/4W, 4300 OHM	1
24	900074	RESISTOR, COMP., CARBON 5%, 1/4W 1500 OHM	1
25	900123	RESISTOR, COMP., CARBON 5%, 1/4W 100K OHM	1
26	900309	RESISTOR, COMP., CARBON 5%, 1/2W, 33K OHM	1
27	66300	RESISTOR, WIREWOUND, 5% 10W, 500 OHM	1
28	63856	TRIAC, 6A, 400V, #SC1410	2
29	66136	POWER TRANSFORMER, P.C. MOUNT, 20 V.C.T., 60 M.A.	1
30	67455	INTEGRATED CIRCUIT, QUAD COMPARATOR, SIGNETICS LM 3394 (SOURCE CONTROL)	1
31	940112	INTEGRATED CIRCUIT, MONO-STABLE, MULTIVIBRATION GATES SN74121N	1
32	66137	COUPLER, PHOTOCELL, LED	2
33	66138	REGULATOR, 3 TERMINAL, POSITIVE, 5V	1
34	65962	SOCKET, RELAY, 11 POSITION	1
35	66280	HOLD-DOWN CUP, RELAY, MIDTEX #139-0029-00	1
36	60971	SCREW, MACH., PAN HD., SLOTTED STL, CD, PL, UNC #6-32X 1/2 lg	2
37	60074	HEX NUT, SST, #6-32	2
38	910143	CAPACITOR, FIXED, .1MF, 4000V (MYLAR)	1

Printed Circuit Board Assembly (GLC-3)

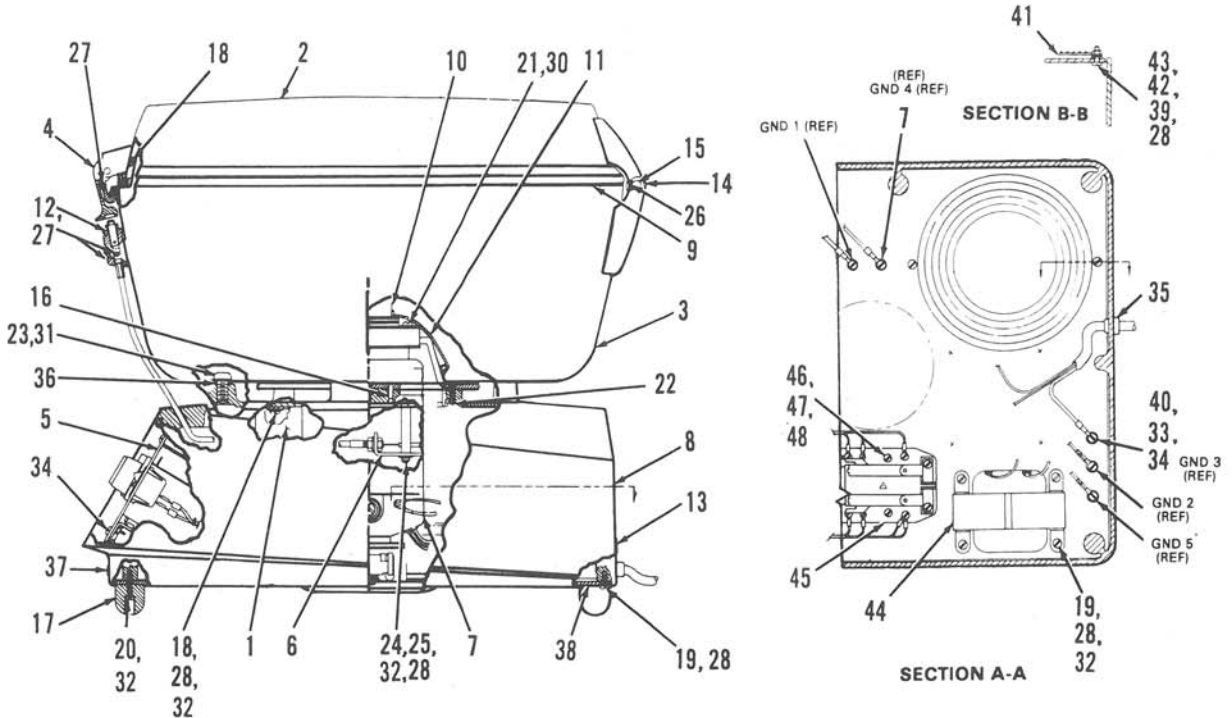

REF.	DWG.NO.	DESCRIPTION	REQ.
1	06107	MICROSWITCH ASSEMBLY	1
2	06188	LID SUB-ASSEMBLY	1
3	06204	GUARD BOND ASSEMBLY	1
4	06190	LATCH SUB-ASSEMBLY	1
5	06294	INSTRUMENT PANEL ASSEMBLY	1
6	06313	LID-LOCK ACTUATOR	1
7	06369	MOTOR ASSEMBLY	1
8	06205	BASE	1
9	06198	COVER SEAL	1
10	06347	GYRO DRIVE ASSEMBLY	1
11	06199	GYRO SEAL	1
12	06143	LOCK BOOT	1
13	06292	NAMEPLATE	1
14	06142	TORSION SPRING	1
15	06186	HINGE PIN	1
16	60237	SPRING PIN, 1/8 x 1/7	1
17	66225	BUMPER, RUBBER, ATLANTIC INDIA #23	4
18	61818	MACHINE SCREW, PAN HD. 8-32 x 5/16	5
19	62358	MACHINE SCREW, PAN HD. 8-32 x 3/8	7
20	64695	MACHINE SCREW, PAN HD. 8-32 x 7/8	4
21	60321	MACHINE SCREW, RD. HD. 10-24 x 1/2	3
22	62246	MACHINE SCREW, FLAT HD. 10-24 x 3/4	4
23	62694	CAP SCREW, BUTTON HD. SOC. 3/8-24 x 3/4	4

REF.	DWG.NO.	DESCRIPTION	REQ.
24	66179	SPACER, ROUND 19/16 LG.	2
25	60152	CAP SCREW, FILL, 8-32 x 17/8	2
26	60173	SET SCREW, CUT POINT, 6-32 x 1/8	1
27	60177	SET SCREW, CUT POINT, 8-32 x 3/16	3
28	60172	LOCKWASHER, SPLIT #8	17
29	61506	LOCKWASHER, INTERNAL, #8	3
30	63346	LOCKWASHER, SPLIT, #10	3
31	61889	NYLON WASHER	4
32	60171	FLAT WASHER, #8	15
33	61711	LOCKWASHER, INTERNAL, #5	5
34	60711	MACHINE SCREW, TRUSS HD., #6-32 x 1/4	9
35	66287	STRAIN RELIEF BUSHING	1
36	62033		4
37	64803	SERIAL NUMBER PLATE	1
38	05341	BASE PLATE	1
39	61652	FLAT WASHER, SST #10	1
40	60522	FLAT WASHER, #6	5
41	66220	FINGER GUARD	1
42	60184	HEX NUT, SST, #8-32	2
43	60964	MACHINE SCREW, PAN HD., SLOTTED, UNC #8-32 x 1/2 Lg.	2
44	66665	FILTER CHOKE, 25MH, SADC	1
45	61002	HEAVY DUTY POWER RELAY, 25A, 12 VAC	1
46	60948	MACHINE SCREW, PAN HD., #8-32 x 5/8 Lg, ST1	2
47	60172	LOCKWASHER, SPLIT, #8	2
48	63085	FLAT WASHER, .1691D	2

General Laboratory Centrifuge (GLC-3)

11/80
GLCS/35-1

Figure 21-1. General Laboratory Centrifuge (GLC-3)
(Sheet 1 of 2)

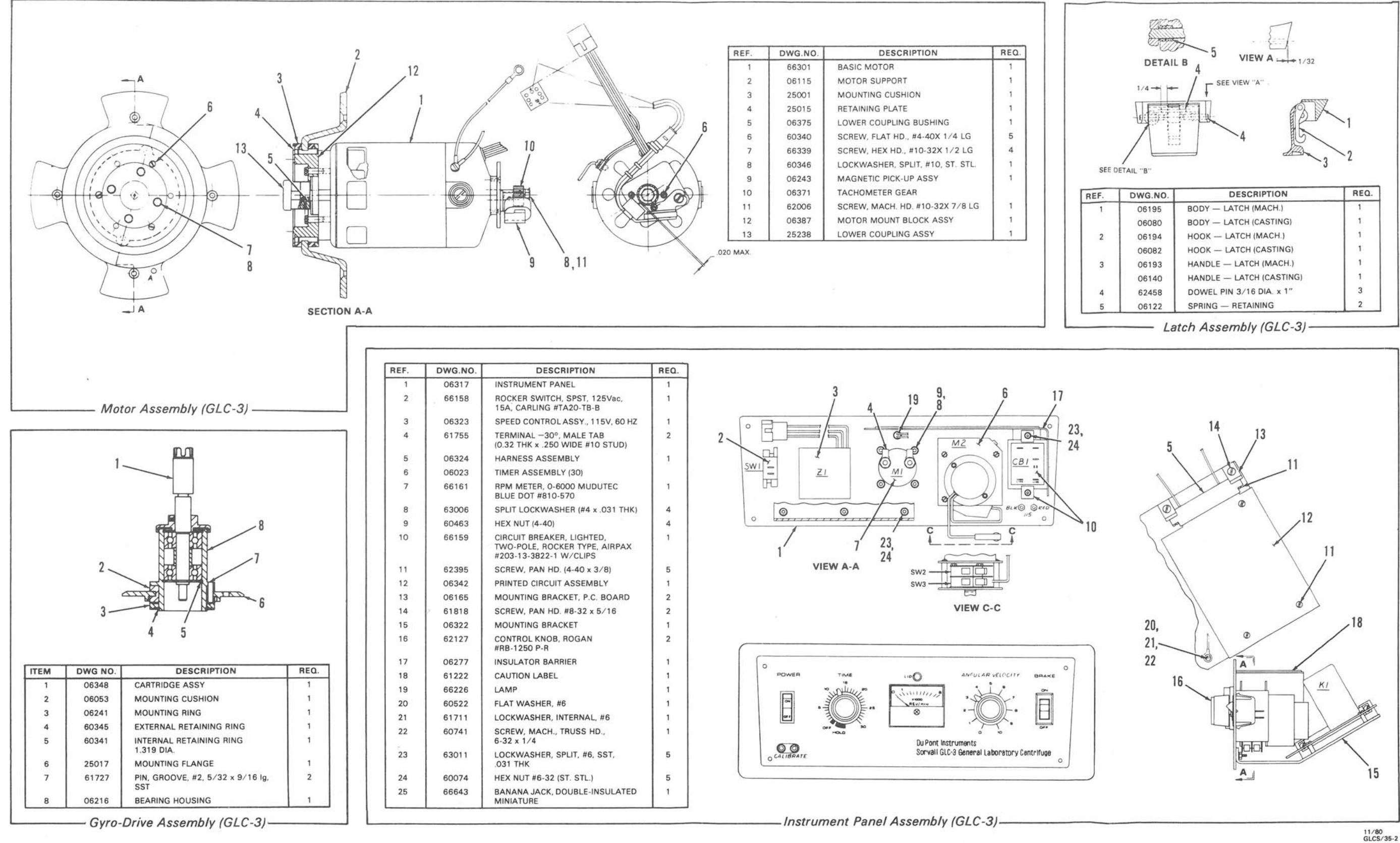


Figure 21-1. General Laboratory Centrifuge (GLC-3)
(Sheet 2 of 2)