



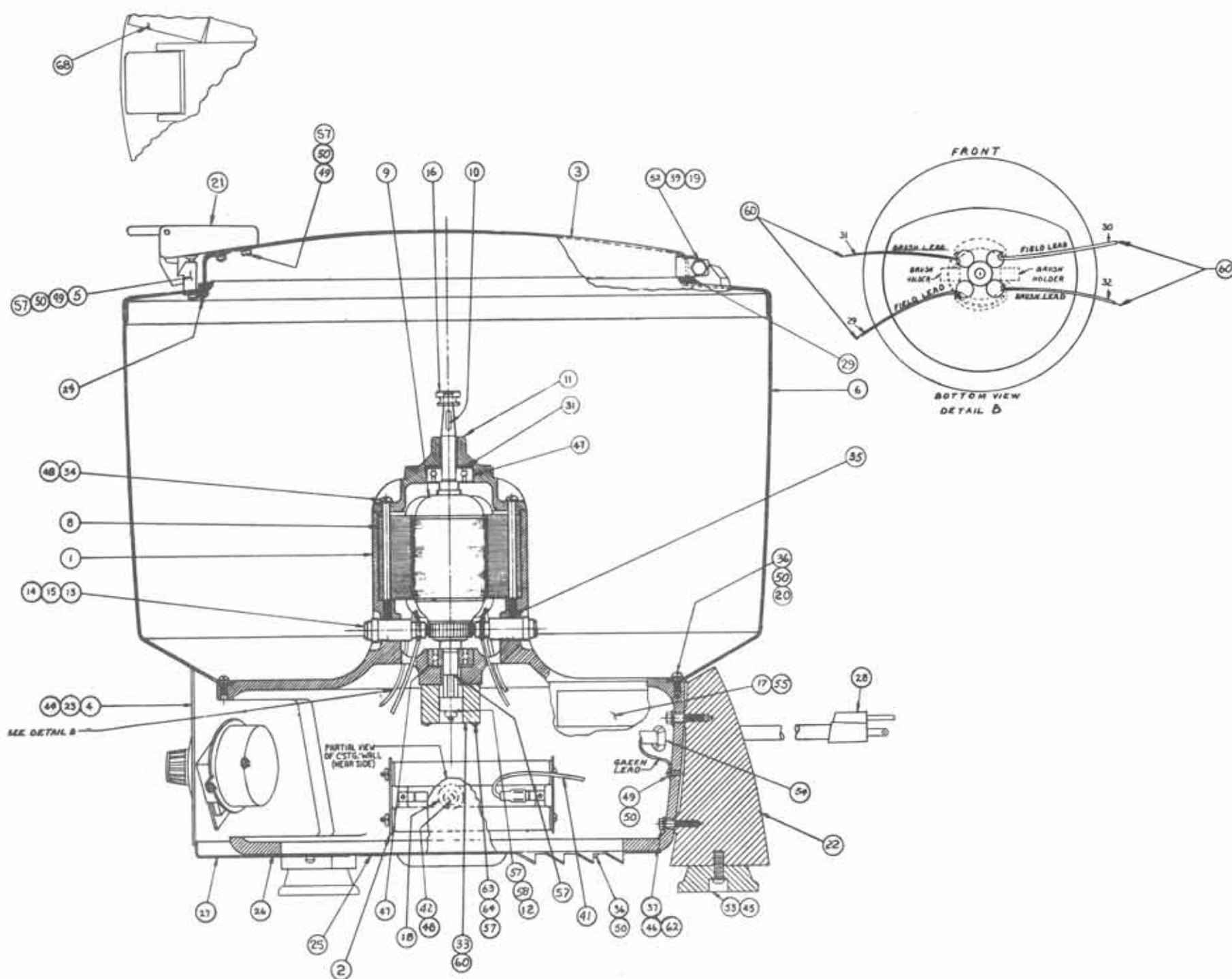

ITEM	DESCRIPTION	PART NO.
1	ASSEMBLY, BASE	42967
2	ASSEMBLY, BRAKE RESISTOR R4	41313
3	ASSEMBLY, GUARD BOWL COVER	46764
4	ASSEMBLY, CONTROL PANEL (SEE FIGURE 5-3)	46753
5	BLOCK CLAMP	46822
6	ASSEMBLY, GUARD BOWL	46765
7	DELETED	----
8	MOTOR FIELD B1	1954
9	MOTOR ARMATURE B1	41101
10	KEY	32385
11	TOP CAP - MACH	42249
12	HOLDER, MAGNET	41100
13	BRUSH HOLDER (TWO) (PT 1)	9037
14	CAP, BRUSH HOLDER (TWO) (PT 2)	9036
15	BRUSH AND SPRING (TWO)	1780
16	NUT, KNURLED	1729
17	PLATE, DATA	30633
18	SPACER, BRAKE RESISTOR	41098
19	LOCKWASHER, STEEL, CAD/PL, 1/4 INT TOOTH (TWO)	COM'L
20	GASKET	42269
21	ASSEMBLY, LATCH	46828
22	LEG, CENTRIFUGE (THREE)	42324A
23	BRACKET (TWO)	32383
24	DELETED	----
25	LABEL, WARNING	45997
26	BOTTOM PLATE	32367
27	GUARD, ELECTRICAL	32390
28	POWER CORD	38805
29	RUBBER CUSHION (FOUR)	1779
30	DELETED	----
31	WASHER, SPRING (WALLACE BARNES #5-10)	41778A
32	DELETED	----
33	TACHOMETER GENERATOR	41461
34	SCREW - RD-HD-SLOT STL CAL PL (TWO) 10-32 x 3.25	COM'L
35	SCREW-SET-SOC-MYLOC-STL CAD PL (TWO) 10-32 x .25	COM'L
36	SCREW - RD-HD-PHIL (TEN) STL CAD PL 10-32 x .38	COM'L
37	SCREW, HEX HD STL CAD PL 1/4-20 x 1.25	COM'L
38	DELETED	----
39	SCREW-SET-SLOT-CONE PT - STL CAD PL 1/4-20 x .62	COM'L
40	DELETED	----
41	WIRING HARNESS, BRAKE RESISTORS	46760
42	SCREW - RD-HD-PHIL STL CAD PL 10-32 x 1.25	COM'L
43	DELETED	----

ITEM	DESCRIPTION	PART NO.
44	SCREW-TRUSS-HD-PHIL STL CAD PL 8-32 x 1.25 (FOUR)	COM'L
45	SCREW-RD-HD-PHIL STL CAD PL 1/4-20 x 1.00 (THREE)	COM'L
46	LOCKWASHER, SPLIT (SIX)	COM'L
47	BEARING (TWO)	9939A
48	LOCKWASHER (FIVE) STL CAD PL #10	COM'L
49	SCREW RD HD PHIL STL CAD PL 8-32 x .32 (SIX)	COM'L
50	LOCKWASHER STL CAD PL #8 (FIFTEEN)	COM'L
51	DELETED	----
52	NUT, ACORN, SMALL, BRASS, CHROME PL 1/4-20 (TWO)	COM'L
53	MOUNTING FOOT (THREE) (ATLANTIC INDIA RUBBER #106)	33340
54	STRAIN RELIEF (HEYCO #SR-6L-1)	COM'L
55	SCREW, DRIVE (TWO) STN STL (#0X.19LG)	COM'L
56	NUT, HEX (TWO)	COM'L
57	LOCTITE (AS REQ.) (COMP #242)	COM'L
58	NUT, HEX STL CAD PL 6-32	COM'L
59	DELETED	----
60	CONTACT, SOCKET (SIX) (AMP #350536-1)	COM'L
61	CLOSED END SPLICE (TWO)	COM'L
62	WASHER, PLAIN STL CAD PL 1/4 (SIX)	COM'L
63	SCREW RD HD SLOT STL CAD PL 4-40 x 1.25 (TWO)	COM'L
64	WASHER, PLAIN (TWO) STL CAD PL #4	COM'L
65	BUTT SPLICE	COM'L
66	DELETED	----
67	ASSEMBLY, JUMPER	46767
68	LABEL, FDA CODE 36980	46290

Figure 5-1 Centrifuge Assembly, 115 V 60 Hz, 2355

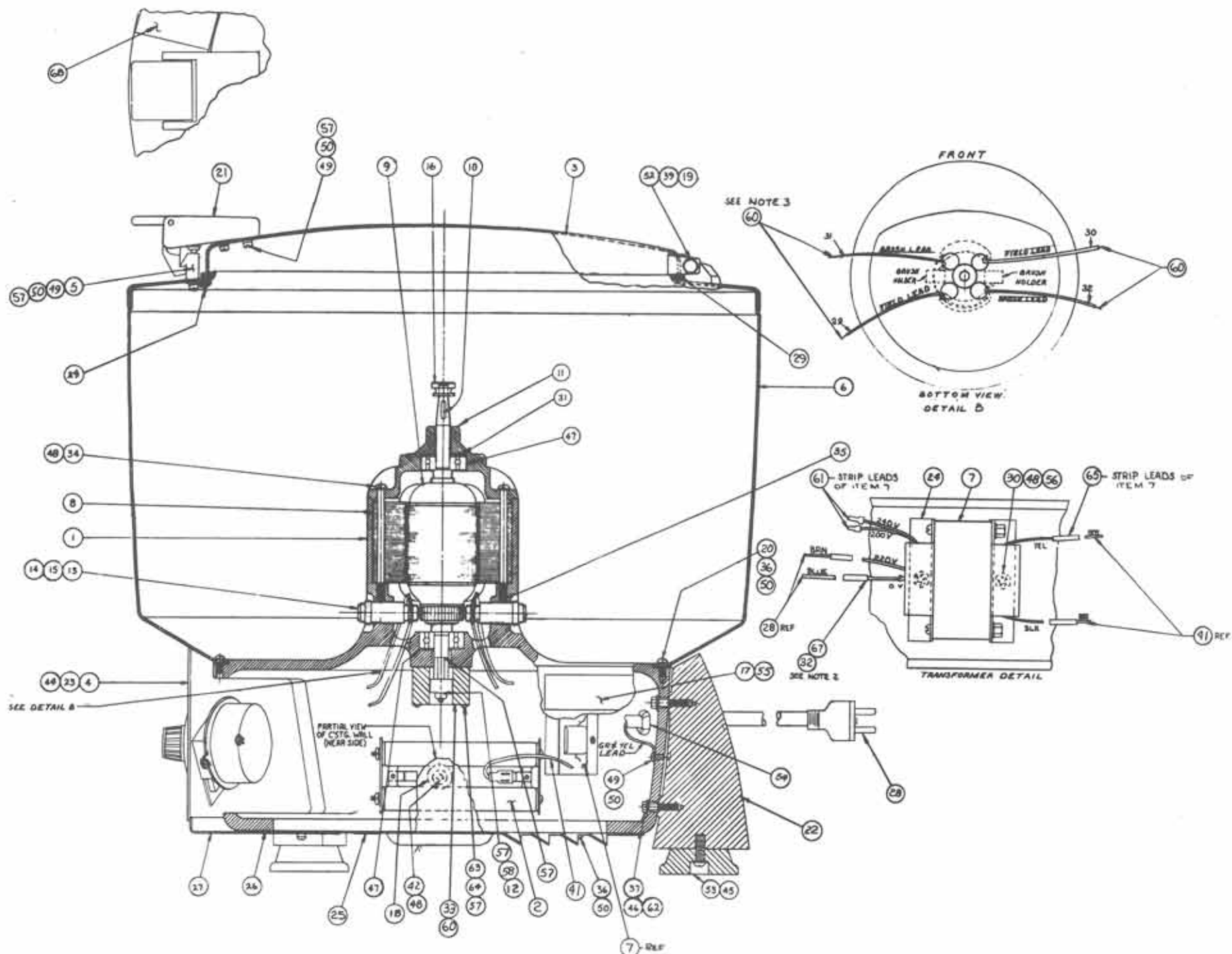

ITEM	DESCRIPTION	PART NO.
1	ASSEMBLY, BASE	41110
2	ASSEMBLY, BRAKE RESISTOR R4	41313
3	ASSEMBLY, GUARD BOWL COVER	46764
4	ASSEMBLY, CONTROL PANEL (SEE FIGURE 5-3)	46756
5	BLOCK CLAMP	46822
6	ASSEMBLY, GUARD BOWL	46765
7	TRANSFORMER	32391
8	MOTOR FIELD B1	1954
9	MOTOR ARMATURE B1	41101
10	KEY	32385
11	TOP CAP - MACH	42249
12	HOLDER, MAGNET	41100
13	BRUSH HOLDER (TWO) (PT 1)	9037
14	CAP, BRUSH HOLDER (TWO) (PT 2)	9036
15	BRUSH AND SPRING (TWO)	1780
16	NUT, KNURLED	1729
17	PLATE, DATA	30633
18	SPACER, BRAKE RESISTOR	41098
19	LOCKWASHER, STEEL, CAD/PL, (TWO)	COM'L
20	GASKET	42269
21	ASSEMBLY, LATCH	46828
22	LEG, CENTRIFUGE (THREE)	42324A
23	BRACKET (TWO)	32383
24	BRACKET, ANGLE (TWO)	32398
25	LABEL, WARNING	45997
26	BOTTOM PLATE	32367
27	GUARD, ELECTRICAL	32390
28	POWER CORD ASSEMBLY EUROPEAN	46759
29	RUBBER CUSHION (FOUR)	1779
30	SCREW, RD HD PHIL (TWO)	COM'L
31	WASHER, SPRING	41778A
32	SOLDER 60/40 (AS REQ.)	COM'L
33	TACHOMETER GENERATOR	41461
34	SCREW - RD-HD-SLOT STL CAD PL (TWO) 10-32 x 3.25	COM'L
35	SCREW-SET-SOC-NYLOC-STL CAD PL (TWO) 10-32 x .25	COM'L
36	SCREW - RD-HD-PHIL (TEN) STL CAD PL 10-32 x .38	COM'L
37	SCREW, HEX HD STL CAD PL 1/4-20 x 1.25	COM'L
38	DELETED	----
39	SCREW, SET-SLOT-CONE PT - STL CAD PL 1/4-20 x .62	COM'L
40	DELETED	----
41	WIRING HARNESS, BRAKE RESISTORS	46760
42	SCREW-RD-HD-PHIL STL CAD PL 10-32 x 1.25	COM'L
43	DELETED	----

ITEM	DESCRIPTION	PART NO.
44	SCREW-TRUSS-HD-PHIL STL CAD PL 8-32 x 1.25 (FOUR)	COM'L
45	SCREW - RD-HD-PHIL STL CAD PL 1/4-20 x 1.00 (THREE)	COM'L
46	LOCKWASHER (SIX) STL CAD PL 1/4	COM'L
47	BEARING (TWO) 4FAFNIR #200 PP	9939A
48	LOCKWASHER (THREE) STL CAD PL #10	COM'L
49	SCREW RD-HD-PHIL STL CAD PL 8-32 x .31 (SIX)	COM'L
50	LOCKWASHER STL CAD PL #8 (FIFTEEN)	COM'L
51	DELETED	----
52	NUT, ACORN, SMALL, BRASS, CHROME PL 1/4-20 (TWO)	COM'L
53	MOUNTING FOOT (THREE)(ATLANTIC INDIA RUBBER #106)	33340
54	STRAIN RELIEF (HEYCO #SR-6L-1)	COM'L
55	SCREW, DRIVE (TWO) STN STL (#OX.19LG)	COM'L
56	DELETED	----
57	LOCTITE (AS REQ.) (COMP #75 OR #59)	COM'L
58	DELETED	----
59	SEALER (AS REQ.) (3M #EC-821)	COM'L
60	CONTACT, SOCKET (SIX) (AMP #350536-1)	COM'L
61	DELETED	----
62	WASHER, PLAIN STL CAD PL 1/4 (SIX)	COM'L
63	SCREW RD-HD-SLOT STL CAD PL 4-40 x 1.25 (TWO)	COM'L
64	WASHER, PLAIN (TWO) STL CAD PL #4	COM'L
65	DELETED	----
66	DELETED	----
67	DELETED	----
68	LABEL, FDA CODE 36980	46290

Figure 5-2 Centrifuge Assembly, 230 V 50 Hz, 2356