



**CENTRIFUGE—EXPLODED VIEW  
CENTRA-MP4R**

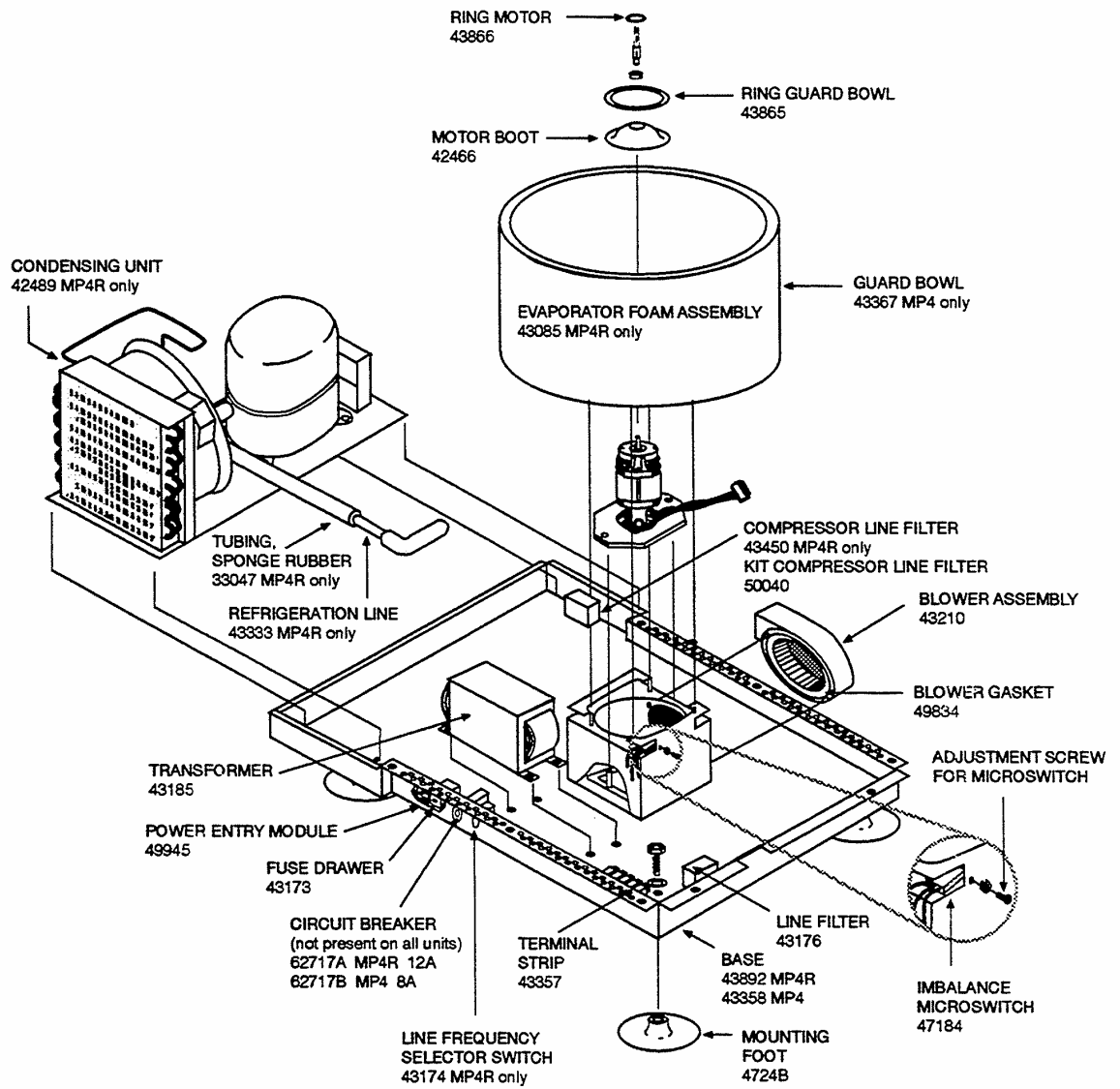


**BASE ASSEMBLY—EXPLODED VIEW  
CENTRA-MP4R**

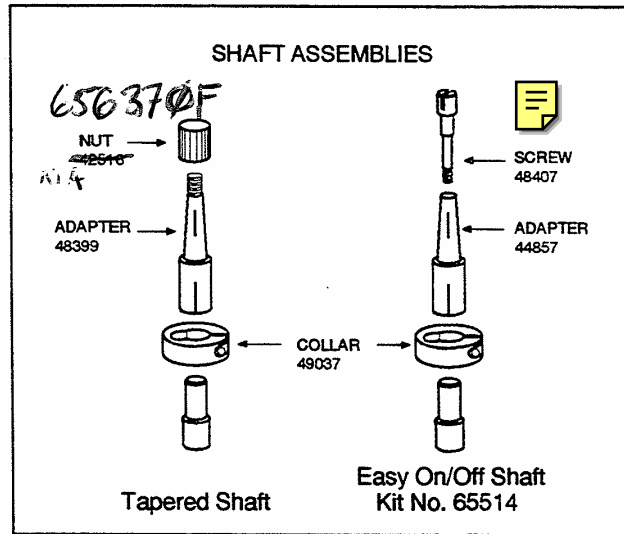

**44990 MOTOR REPLACEMENT KIT (TO REPLACE 43310 WITH 44996)**

Software versions earlier than 2.0 is 43310 Motor  
 Software versions 2.0 or later is 44996 Motor  
 To verify version level see Section 10.4 Special Functions.

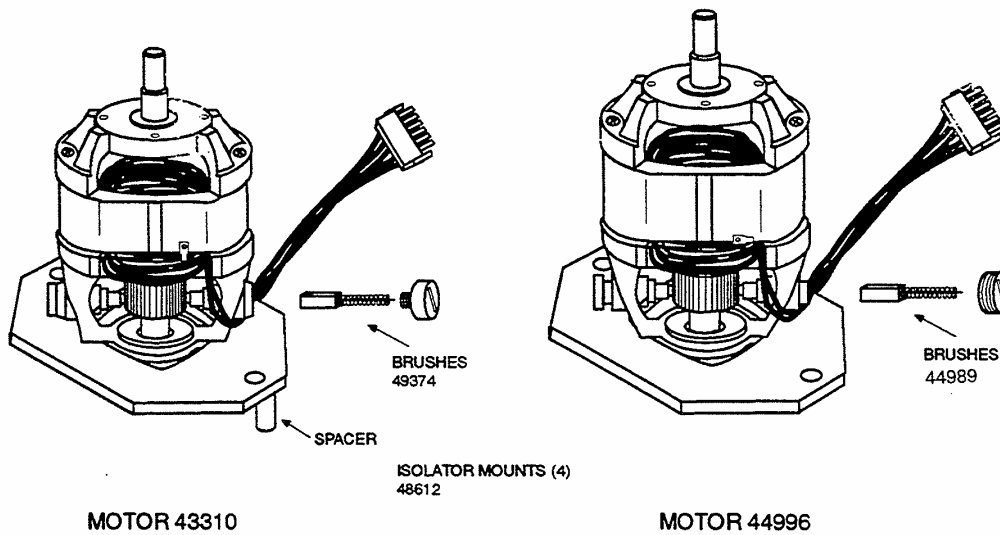

**CENTRA-MP4/R DRIVE**

**\*Assy Includes**

43304        Motor Assy  
    43310        Motor  
    43870        Blower Mount Plate  
    33065        Channel Pawling  
    43125        Drive Frame  
    43135        Cylinder  
    43150        Cover Plate  
    43195        Imbalance Harness  
    43210        Blower Assy  
    43864        Motor Plate Box  
    47184        Microswitch  
    48612        Isolator  
    49834        Blower Box Gasket

43706        Spare Parts Kit  
    43689        Fuse 4A - 240V (4)  
    43707        Fuse 8A - 120V (4)  
    43177        Line Cord  
    43312        Line Cord (International)  
    49374        Brushes  
    65436        MAG. SOCKET

Centra MP4 2437

\*see included parts list

43358	Base Plate	
43210	Blower Assy	
49374	Brushes	
43348	Control Panel (no PCB)	
49847	Control Panel Clip	
45249	Cabinet	
65253	Cover Bumper	E-From 50006
40502	Cover Gasket	
65435	Cover Insert	
43381	Cover Insert Plate	QA CIRCUIT BREAKER 62717B
43890	Cover Panel	
47248	Foot-Rubber	
43174	Frequency Selector	
43689	Fuse 4A - 240V	
43707	Fuse 8A - 120V	
49937 43250	Fuse PCB 2.5A (F1)	
9946	Fuse PCB 6.3A (F2)	
43173	Fuse Drawer	
43367	Guard Bowl	
43641	Harness Blower/PCB	
43195	Harness Imbalance	
43253	Harness Latch/PCB	
43254	Harness Motor/PCB	
43255	Harness P.E.M.	
43251	Harness Power/PCB	
49825A	Hinge	
47184	Imbalance Switch	N/S 50115
43181A	Latch	
43181B	Latch Striker	
43177	Line Cord	
43176	Line Filter	
43310	Motor	
43304	Motor Assy*	
42466	Motor Boot	
44586	P.C. Board	
9555	Plug Bypass Hole	
43175	Power Entry Module	
43261	Rear Access Panel	
50242-43866	Ring-Motor	BULK PLASTIC
43865	Ring-Guard Bowl	
42516	Rotor Locking Nut	
48399	Shaft Adapter-Tapered	E-Z ON SHAFT ADAPTER KIT 65514
49037	Shaft Adapter Collar Clamp	
43706	Spare Parts Kit*	SCREW KIT 50308
43185	Transformer	
48110	Viewport	