

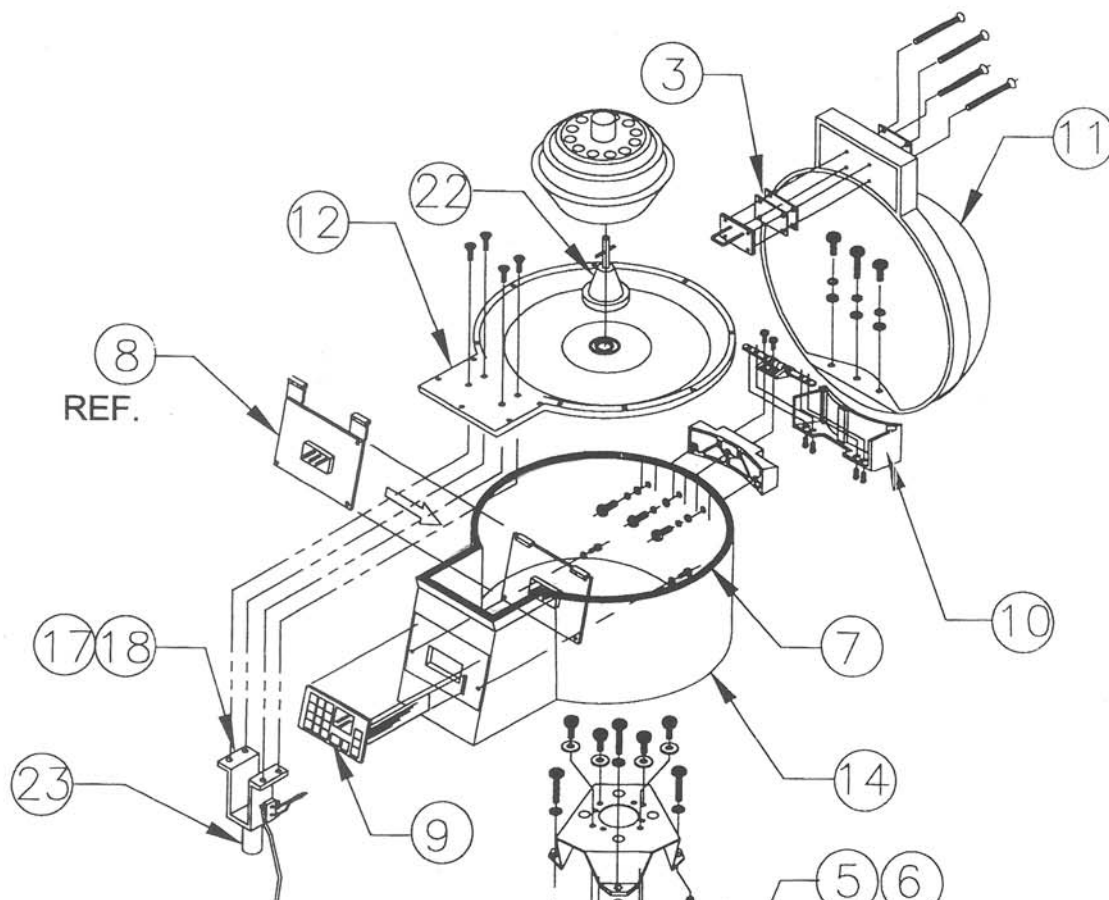


SEROFUGE R/P LIST

ITEM	PART DESCRIPTION	420351	420352	420354
1	FUSE, 115 V	421338	421338	N/A
2	FUSE, 230V	N/A	N/A	421339
3	LATCH HOOK KIT	421340	421340	421340
4	PEM ADAPTER MODULE,115V	421341	421341	421341
5	POWER ENTRY MODULE,115V	421342	421342	N/A
6	POWER ENTRY MODULE,230V	N/A	N/A	421355
7	LID SEAL	421343	421343	421343
8	PCB ASSEMBLY KIT	421344	421344	421344
9	KEYPAD	421345	421354	421345
10	HINGE SET KIT	422953	422953	422953
11	LID	421347	421347	421347
12	ROTOR BOWL	421348	421348	421348
13	BASE	421349	421349	421349
14	SHELL	421350	421350	421350
15	MOTOR KIT, 115V	421351	421351	N/A
16	MOTOR KIT, 230V	N/A	N/A	421353
17	LATCH ASSEMBLY KIT, 115V	421359	421359	N/A
18	LATCH ASSEMBLY KIT, 230V	N/A	N/A	421356
19	12 PLACE ROTOR	420545	420545	420545
20	8 PLACE ROTOR	420224	420224	420224
21	6 PLACE ROTOR	420526	420526	420526
22	SPINDLE ASSY	421367	421367	421367
23	SOLENOID	429599	429599	429600

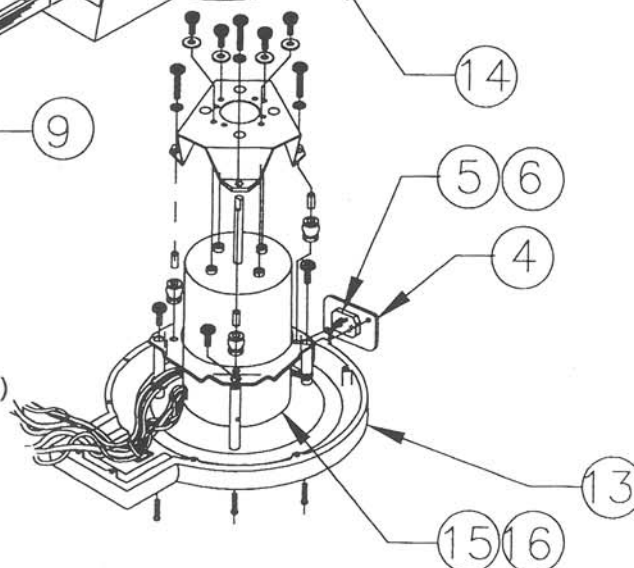

WIRE LIST			
PIN NO.	WIRE COLOR	PIN NO.	WIRE COLOR
1	BLK	8	BLK
2	WHT	9	BLK
3	GRN	10	WHT
4	BLK	11	RED
5	BLK	12	BLK
6	YEL/GRN	13	BLK
7	BLK	14	BLK

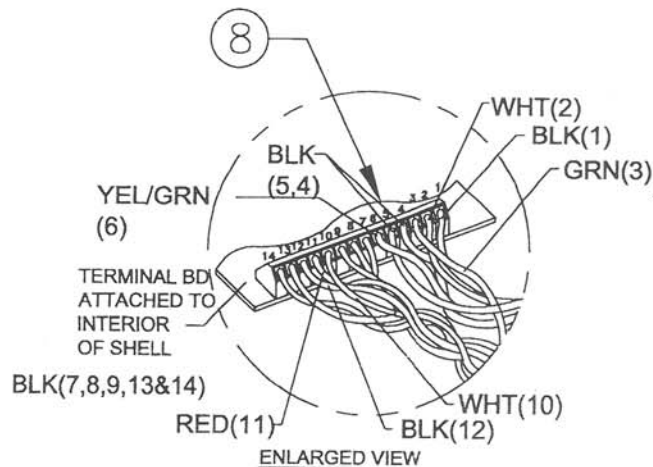

TITLE: SEROFUGE I.P.B
R/P GUIDE

AW-0350-000-078

REV. B