



6 SPARE PARTS LIST

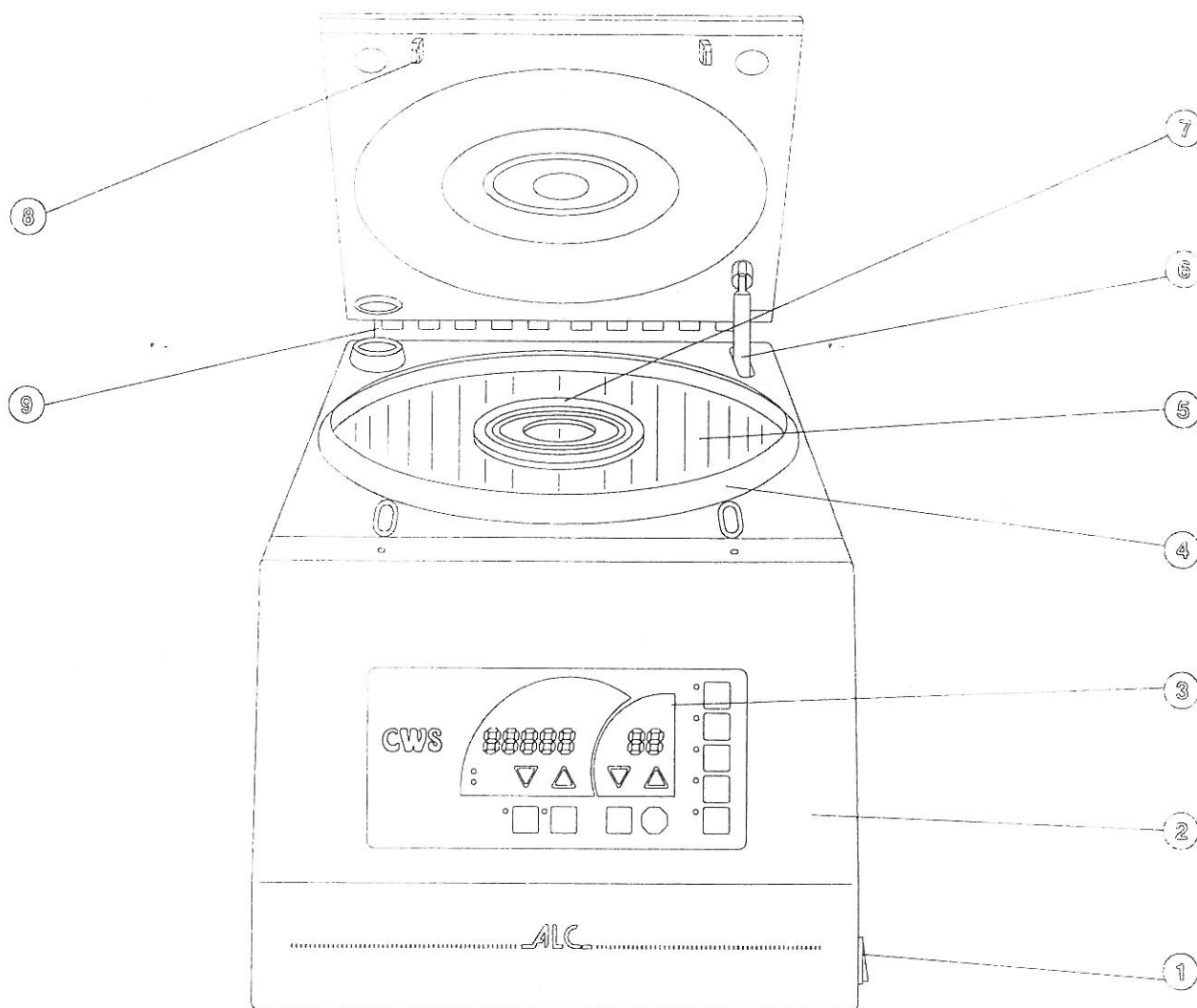


GENERAL VIEW

ITEM	Catalog no.	Description	Qty.
1	89000838	Power switch	1
2	34033002	Front Panel	1
3	34370015	Keyboard, overlay	1
4	89000943	Bowl seal, lid	1
5	89002039	Centrifuge chamber	1
6	89000840	Gas spring	1
7	89000839	Bowl seal chamber, motor	1
8	85240195	Lid latch catch	2
9	11200717	Lid hinge	1
	85280515	Main P.C.B.	1
	11202402	Display P.C.B.	1

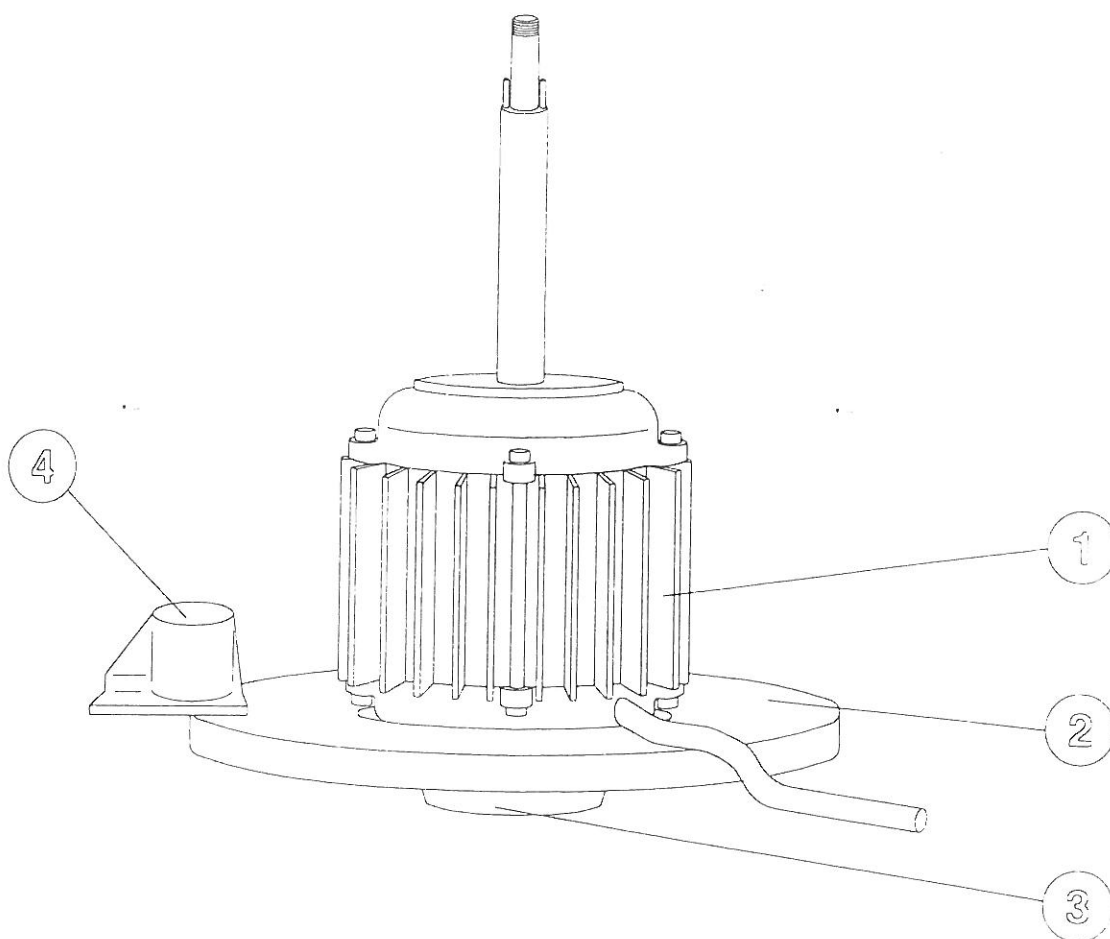


MOTOR

Item	Catalog No.	Description	Qty.
1	38500012	Motor	1
2	11200508	Counter Weight	1
3	89000853	Tachometer Sensor Assy. + Shock Mount	1
4	89000843	Load Imbalance Sensor	1

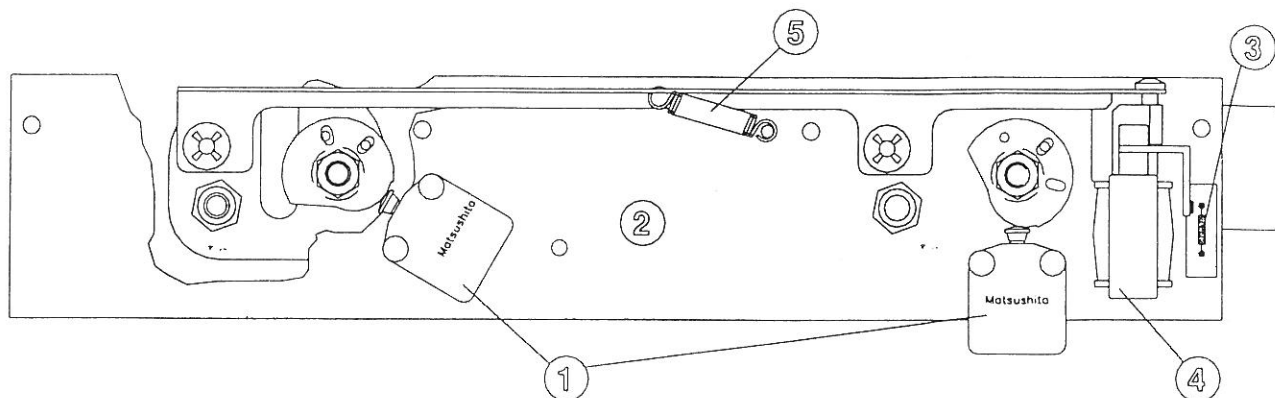

LOCK VIEW

Item	Catalog No.	Description	Qty.
1	89000842	Lid microswitch	2
2	11202913	Lock Assembly	1
3	11250006	Solenoid switch	1
4	11250002	Lid solenoid kit	1
5	89000846	Spring	1

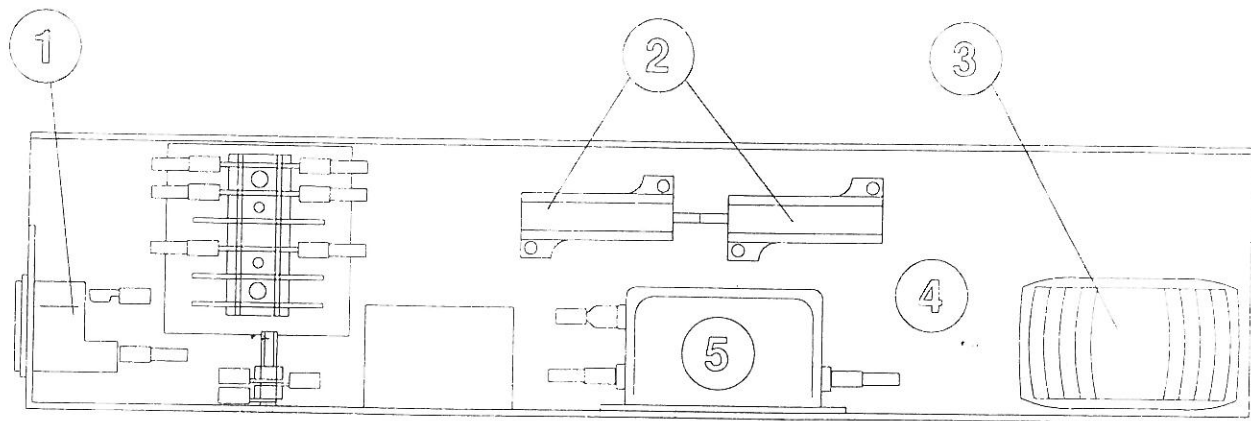

POWER RACK

Item	Catalog No.	Description	Qty.
1	86000775	Mains receptacle	1
2	86003247	Braking resistor	2
3	86003242	Transformer, 230V	1
3	86003243	Transformer, 120V	1
4	11202935	Power supply rack, 230V	1
4	11202936	Power supply rack, 120V	1
5	86003248	Line filter, 230V	1
5	86003256	Line filter, 120V	1