


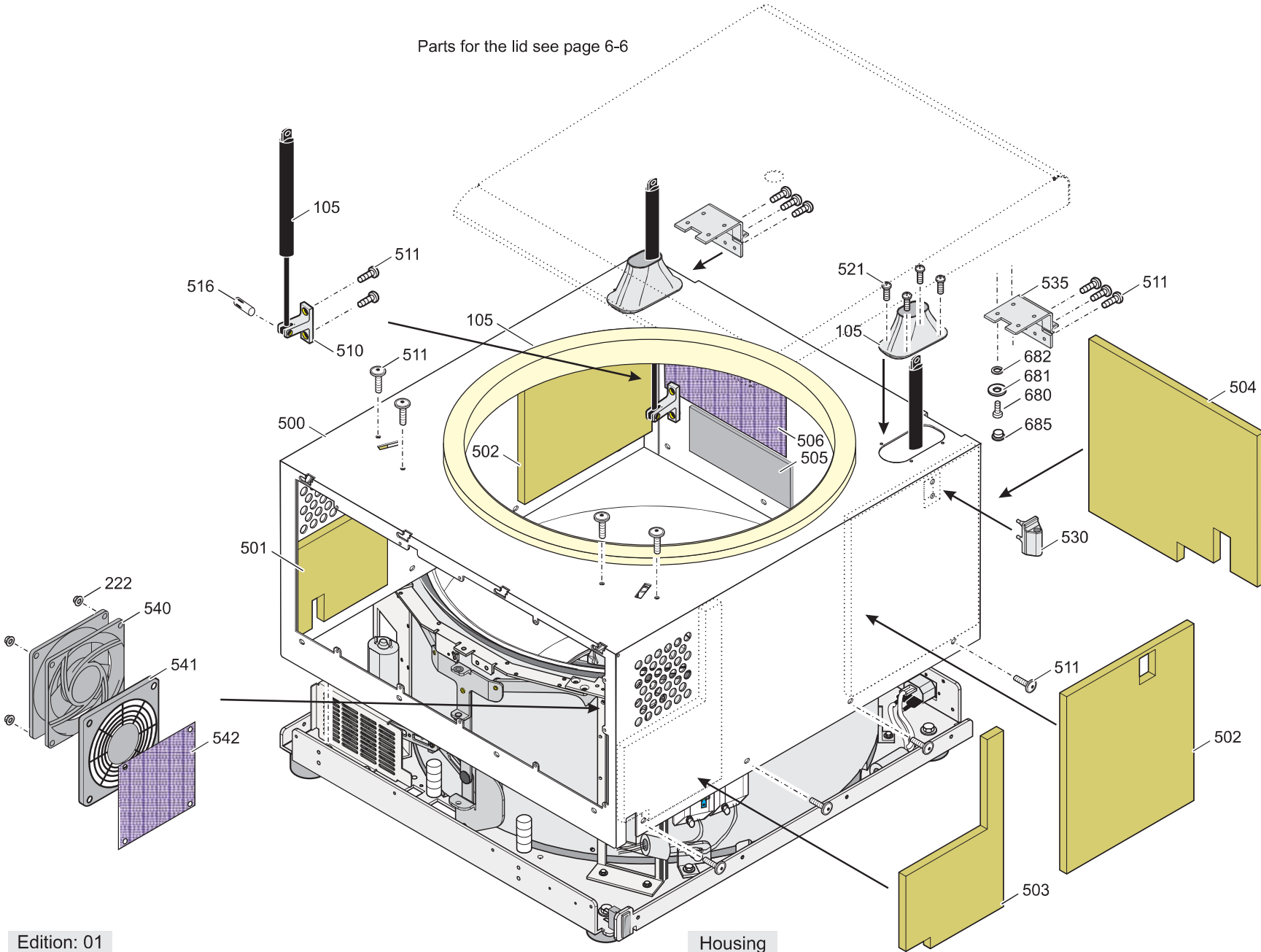



Parts for the lid see page 6-6

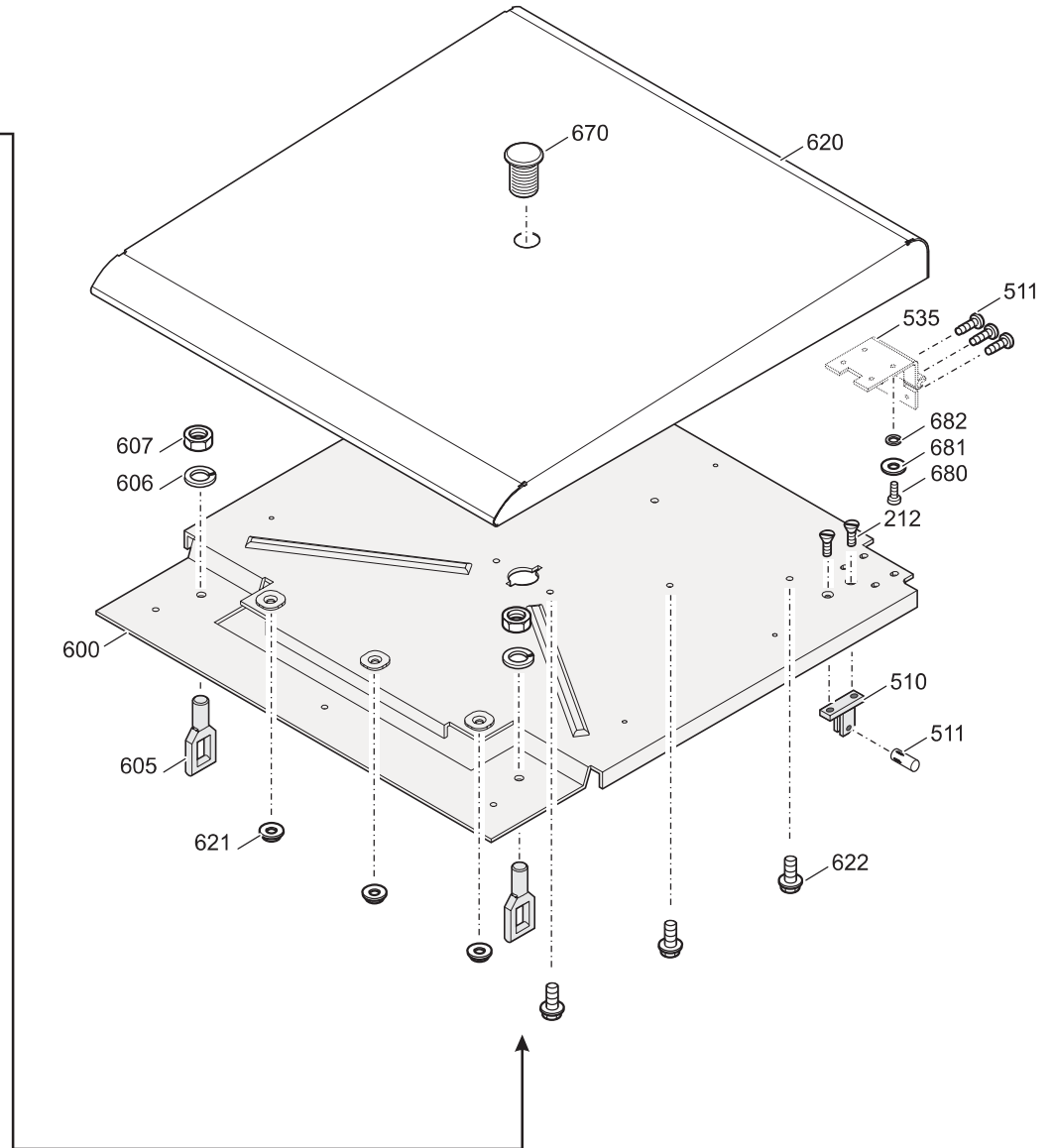

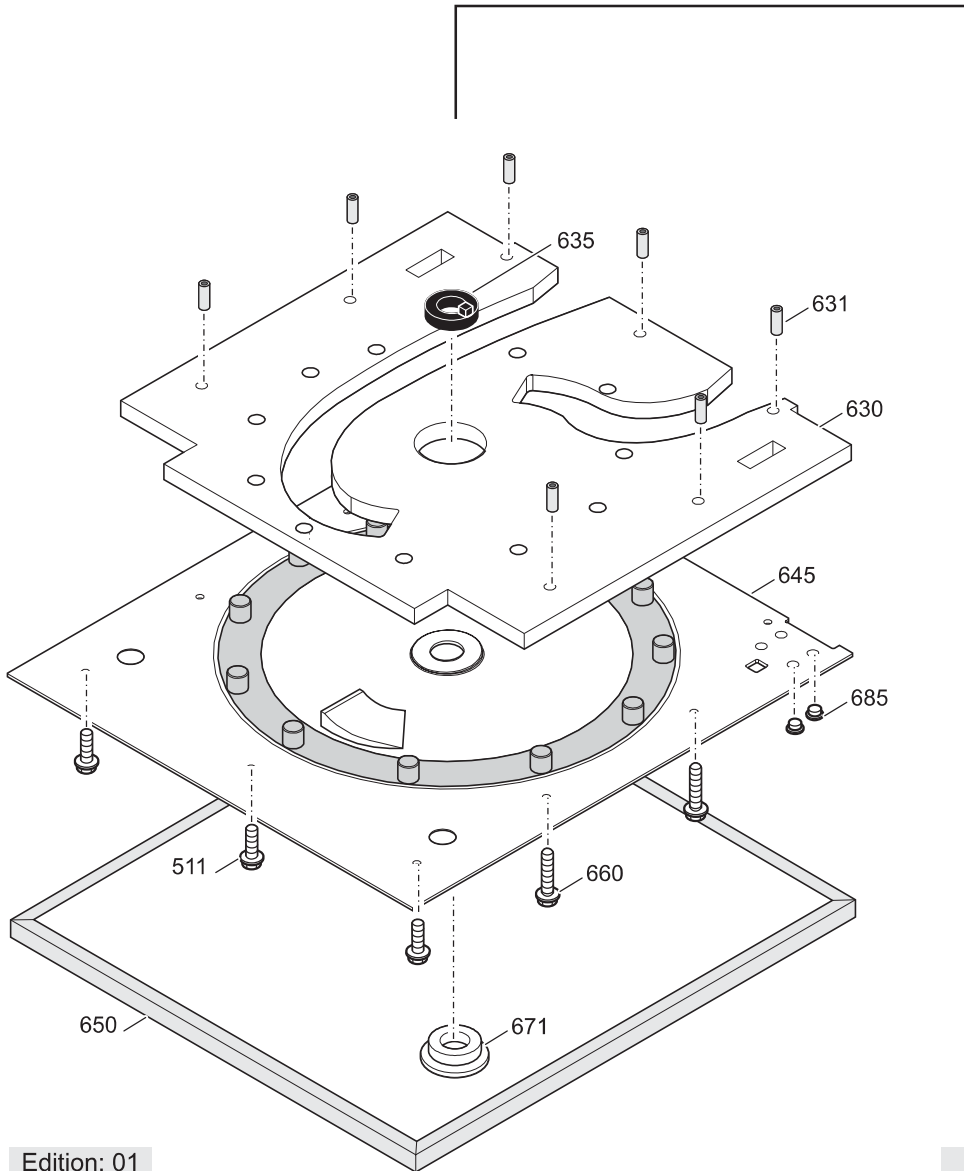

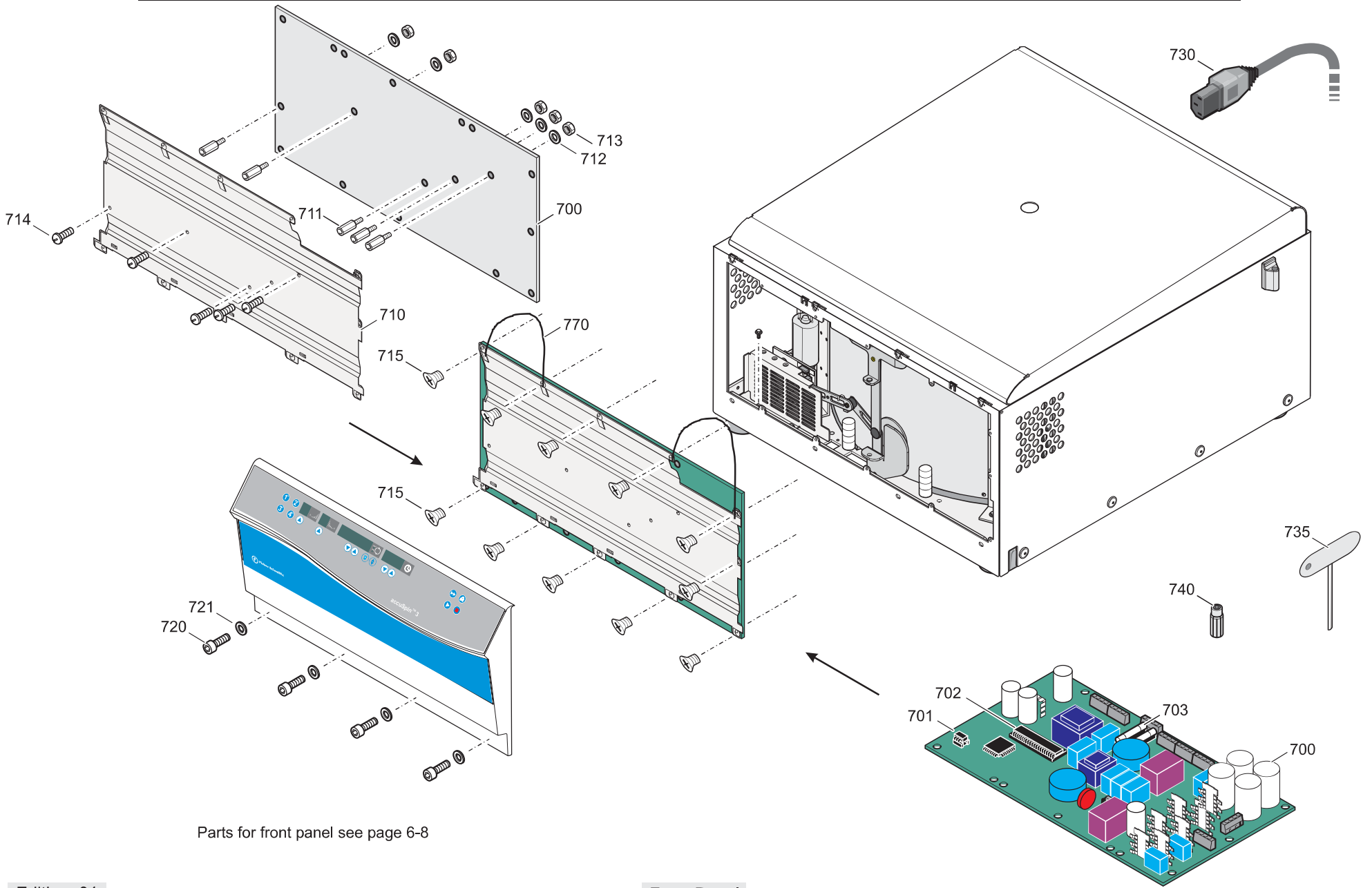

Parts for front panel see page 6-8

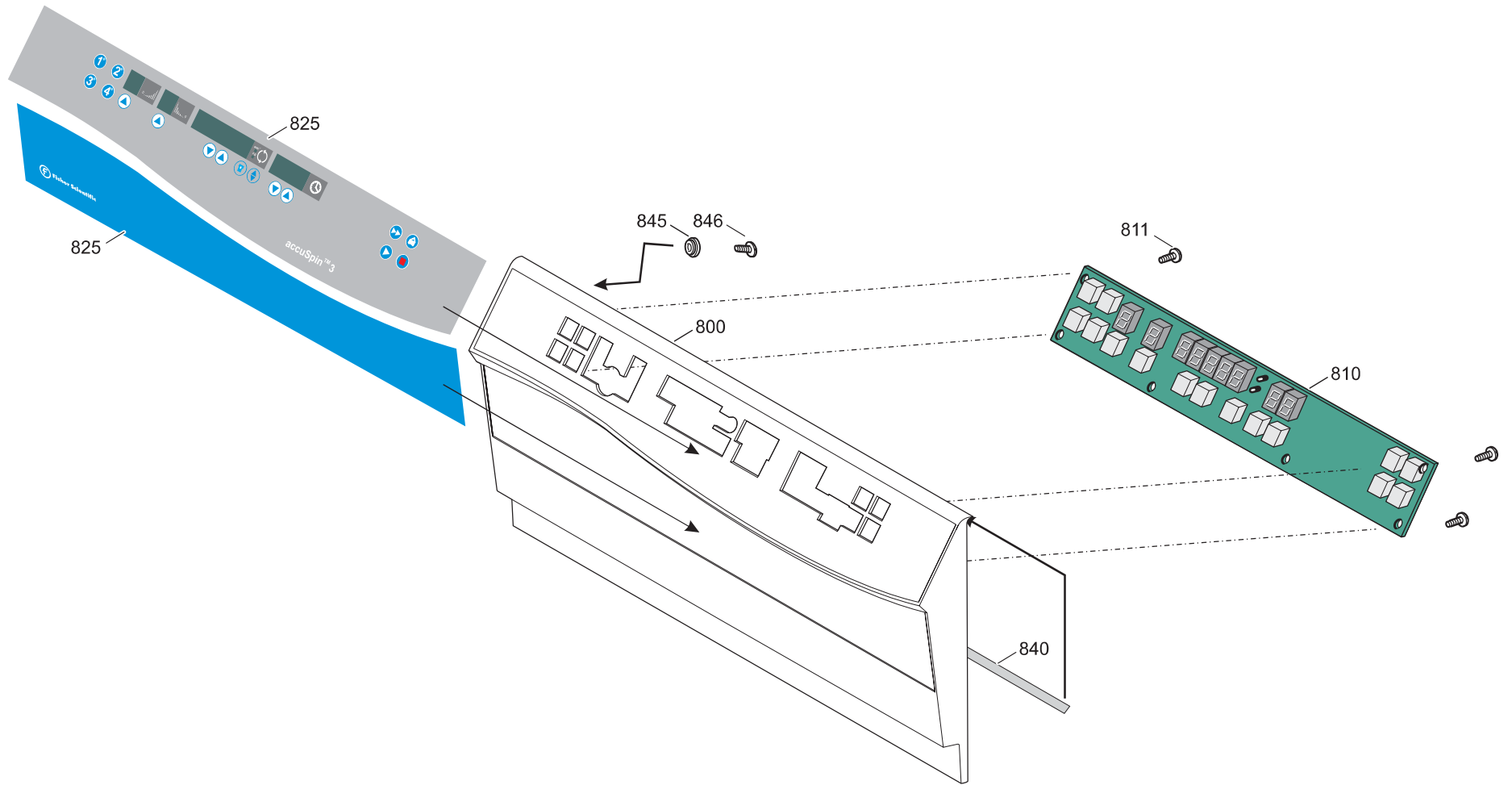

75004391 _01

FISHER ACCUSPIN 3 120 V / 60 HZ

from Serial-No.

to Serial-No.

Index	Partno.	Text
00100	70056750	Chassis
00101	20460089	SCREW M5X5
00102	20057535	Motorchassis
00105	70902661	Parts f maintenance Multifuge/Legend air
00106	20420167	NUT
00109	20510292	SCREW
00110	20430316	Bolt M 12x30
00111	20480247	SECURING RING
00120	20300084	FOOT
00121	20480254	SECURING WASHER
00122	20430041	SCREW M 8 X 12
00125	70057541	Chassis cover
00126	20480244	LOCK WASHER
00127	20460094	SCREW M 4 X 8
00130	70057534	Motorchassis
00135	20210390	MOTOR 125/45 3PH
00137	20151180	Rotorrecognition board JST
00138	20510350	Bolt
00140	20151214	PCB Speed and Imbalance
00141	20460097	Z-SCREW M3 X 8
00142	20480242	SECURING WASHER
00150	20180608	RECEPTACLE
00151	20440023	SCREW M3 X 8
00152	20460090	SCREW M4X12
00153	20420076	HEXAGONAL NUT M4
00155	70060308	Sheet
00160	20180138	CONNECTOR
00161	20480108	FAN WASHER A 4,3
00180	70060313	Bracket
00185	20180683	SWITCH 16A/250V
00200	20290658	SEAL 12 x 10
00210	20170330	FERROXCUBE
00222	20420165	NUT
00225	20210401	AXIAL FAN TYP 4606TA 115V 50/60HZ
00250	20056764	Safety chamber
00251	20510353	Screw
00255	20160733	RESISTOR 120 OHM
00256	20160734	RESISTOR 120 OHM w Thermoswitch

75004391 - 01

FISHER ACCUSPIN 3 120 V / 60 HZ

from Serial-No.

to Serial-No.

Index	Partno.	Text
00257	20510352	Screw
00258	20170329	Ferroxcube ring
00270	70904553	bowl
00271	20420026	NUT
00310	20290653	Seal
00325	20056761	MOTORCOVER
00340	70057552	Metall sheet
00400	20210391	MOTOR W GEAR
00405	20056715	Holder
00406	20460163	Screw
00410	20150186	PCB SENSOR
00415	70904339	COUPLING
00416	20510288	Set screw
00417	20440285	SCREW
00420	70056790	Sheet, left
00425	70056789	Sheet, right
00430	20056737	Slideplate
00431	20460092	ALLEN SCREW M4 X 6
00432	20290586	SEAL
00440	20056738	Roller
00441	20056739	Bolt 4
00442	20420107	NUT M 6 LEFT
00450	20056792	HOLDER LATCH PROTECTION
00451	20350242	Spring
00452	20056791	LATCH PROTECTION
00453	20460202	Screw Ejot
00454	20057932	Insulation
00460	20180677	SWITCH
00461	20460157	Bolt M 2,3
00465	70056718	Bracket, Latch-
00470	20180654	SWITCH
00471	20057257	Holder sheet
00475	70904336	LATCH ASSY
00476	20350241	Clip
00480	20056707	Bar
00481	20420070	NUT M 6
00500	70056751	Housing
00501	20057390	Insulation, left front

75004391 _01

FISHER ACCUSPIN 3 120 V / 60 HZ

from Serial-No.

to Serial-No.

Index	Partno.	Text
00502	20057391	Insulation, back
00503	20057389	Insulation 210x245right front
00504	20057392	Insulation back right
00505	20057393	Insulation back left
00506	20057482	GRILL PLASTIC
00510	20056650	HOLDER
00511	20510285	SCREW M5 X 12
00516	20490287	PIN
00521	20510294	SCREW M3 X 10
00530	20057553	Holder, Tool-
00535	20056736	HINGE
00540	20210402	AXIAL FAN TYP 3900M 115V 50/60HZ
00541	20210406	Grill
00542	20057394	GRILL
00600	70056741	Insert
00605	20057217	HOOK
00606	20480236	SECURING WASHER
00607	20420121	NUT M 10 X 1,25
00620	70902580	Lid
00621	20420201	Nut
00622	20510352	Screw
00630	20056786	Insert
00631	20056762	Spacer
00635	20057563	Spacer
00645	20056785	Lid lower part
00650	20057833	Seal
00660	20510360	Screw M5
00670	20056732	Window
00671	20056625	WINDOW, LOWER PART
00680	20510021	SCREW M 5 X 12
00681	20480200	WASHER
00682	20480167	FAN WASHER A 5,1
00683	20310504	Cover
00700	20151184	PCB mainboard
00701	70904591	NV RAM accuspin 3
00702	70000528	PROCESSOR MAIN 0528V09
00703	20230172	FUSE 15 A
00710	70060305	Cover

75004391 _01**FISHER ACCUSPIN 3 120 V / 60 HZ**

from Serial-No.

to Serial-No.

Index	Partno.	Text
00711	20500200	Spacer
00712	20480300	Washer M4
00713	20420200	Nut M4
00714	20470263	Screw
00715	20440284	SCREW M 4 X 8
00720	20510293	Screw
00721	20480130	WASHER 4,3
00730	20190300	Cordset NEMA 5-15P 15A/125V 3x14AWG
00735	20360073	SCREW DRIVER
00740	20904430	COLLET CHUCK
00770	20310508	String w washer
00800	20058192	Frontpanel
00810	20151188	PCB Display
00825	20058205	OVERLAY Multi 3.0 120V/60Hz FISHER
00840	20290664	seal
00845	20057560	Spacer
00846	20460161	SCREW