

EBA 280

TABLETOP CENTRIFUGE

SEPARATING OURSELVES 90°
FROM THE COMPETITION

Hettich
CENTRIFUGES

Compact & Powerful Clinical Tabletop Centrifuge

High-performance with a 90° swing-out rotor for flat separation of all standard blood, urine and pediatric tubes. Achieve consistent platelet counts less than 10,000 platelets/ μ L in 3 minutes or less.

Ergonomic Quick-Lift Rotor System

Provides effortless sample management and installation of centrifuge rotor. The rotor is self-standing allowing you to load samples at the benchtop level. Quickly and easily place the auto-locking rotor onto motor shaft for safe centrifugation. Upon completion, simply lift the rotor out of centrifuge for convenient unloading of samples. No tools or physical stress required.

Memory Control System

Up to 10 pre-programmable settings for easy recall of application parameters. This ensures repeatability and limits programming errors.

Run-Silent™ Operation

Excessive noise in the laboratory can lead to distractions, stress, anxiety, high blood pressure and chronic fatigue. Our centrifuges are as quiet as a normal conversation and will not increase the noise levels within your lab.

Easy-Lift™

Our Easy-Lift Lid, featuring stay-in-place positioning and lid drop protection, is ergonomically designed to limit repetitive stress injuries, spillage and accidental sample disruption when loading and unloading your centrifuge.

Safety-Lock™

Our Safety-Lock Lids are self-closing, auto-locking and come with either single or dual steel lid locks. This ensures a safe, tight closure of the centrifuge chamber during operation. The centrifuge will not operate if lid is not secure. For added safety, our centrifuges feature **all-metal containment** and adhere to all national and international safety standards and certifications.

Five (5) Year Warranty


FDA registered centrifuge manufacturer



MAX CAPACITY:
6 x 50 mL



MAX RPM/RCF:
6,000/5,071

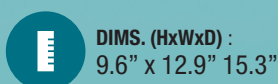

DIMS. (HxWxD) :
9.6" x 12.9" x 15.3"


→ www.hettweb.com/eba280

HETTICH EBA 280 CENTRIFUGE

APPLICATION PACKAGES


All Hettich centrifuge packages come complete with everything you need to perform your application – centrifuge, rotor, buckets, adapters, inserts, bio-containment lids (optional where applicable), five (5) year warranty and free support.

EBA 280 BLOOD TUBE PACKAGE - 6

	Tube Qty	Volume (mL)	Size (mm)	EBA 280	EBA 280S	EBA 280	EBA 280S
				RPM		G-FORCE	
	6	1.6-5	13 x 75	4,700	6,000	2,618	4,266
	6	4-7	13 x 100	4,700	6,000	3,112	5,071
	6	4-7	16 x 75	4,700	6,000	2,865	4,669
	6	8.5-10	16 x 100	4,700	6,000	3,112	5,071


PKG. NO. EBA 280
280BLOOD6
PKG. NO. EBA 280S
280SBL00D6

EBA 280 BLOOD TUBE PACKAGE - 8

	Tube Qty	Volume (mL)	Size (mm)	RPM	G-FORCE
				8	1.6-5
	8	4-7	16 x 75	5,000	2,991


PKG. NO.
280BLOOD8

EBA 280 BLOOD TUBE PACKAGE - 12

	Tube Qty	Volume (mL)	Size (mm)	RPM	G-FORCE
				12	1.6-5
	12	4-7	13 x 100	6,000	4,146
	12	4-10	16 x 75/ 100	6,000	4,146


PKG. NO.
280BLOOD12

EBA 280 PEDIATRIC TUBE PACKAGE - 6

	Tube Qty	Volume (mL)	Size (mm)	EBA 280	EBA 280S	EBA 280	EBA 280S
				RPM		G-FORCE	
	6	0.5	10.7 x 36	4,700	6,000	1,877	3,059


PKG. NO. EBA 280
280BLOOD6-PED
PKG. NO. EBA 280S
280SBL00D6-PED

EBA 280 SEROLOGY PACKAGE - 12

	Tube Qty	Volume (mL)	Size (mm)	RPM	G-FORCE
				12	5

PKG. NO.
280SEROLOGY

EBA 280 SEROLOGY DECANTING PACKAGE - 12

	Tube Qty	Volume (mL)	Size (mm)	RPM	G-FORCE
				12	5


PKG. NO.
280SEROLOGY-D

EBA 280 URINALYSIS PACKAGE - 6

	Tube Qty	Volume (mL)	Size (mm)	RPM	G-FORCE
				6	12 (KOVA)

PKG. NO.
280URINALYSIS6

EBA 280 CUSTOM PACKAGE

	Tube Qty	Volume (mL)	Size (mm)	RPM	G-FORCE
				<p>Additional packages and configurations are available for this device, please inquire. We offer a complete centrifuge product range from small microliter centrifuges, compact tabletop centrifuges, medium to large benchtop centrifuges, high capacity benchtop and floor standing centrifuges for general purpose and specialty applications, inquire for more details.</p>	

PKG. NO.

→ Learn more at: www.hettweb.com/eba280



Discover more solutions online at: www.hettweb.com/centrifuges
Hettich Lab Technology | North America | Toll Free: 1 (866) 370 4388 Email: info@hettweb.com

